

Alt-R™ CRISPR-Cas9 sgRNAs

Chemically synthesized and modified single guide RNAs for outstanding CRISPR performance and quality



**sgRNAs in days,
not weeks,** with fast
synthesis time
(3–5 business days*)



Guaranteed function
with predesigned
sgRNAs



**Custom features to
meet your needs,**
such as a variety of
deliverable sizes,
chemical modifications,
and purification



**Trusted quality and
manufacturing,**
delivering optimized
synthesis and purification
to mitigate oligo cross
contamination risk

* 3–5 business days for most standard requests. Custom requests may require additional manufacturing time.

Alt-R CRISPR-Cas9 single guide RNAs (sgRNAs) comprise both crRNA and tracrRNA sequences within a single molecule. High editing levels are observed at >95% of sites in Jurkat cells (**Figure 1**). Alt-R Cas9 sgRNAs are ideal for challenging conditions such as high nuclease environments or when co-delivered with Cas9 mRNA. They contain chemical modifications that provide increased stability, potency, and resistance against nuclease activity (**Figure 2**) [1].

Customizable sgRNAs to fit every project and every budget

Available in a wide range of deliverable yields, Alt-R Cas9 sgRNAs can be customized to suit small and large experiments. They are available in tube or plate format in a variety of scales from 2 nmol and up. Further options for custom chemical modifications, additional purification, and custom formulation provide flexibility to meet your experimental needs.


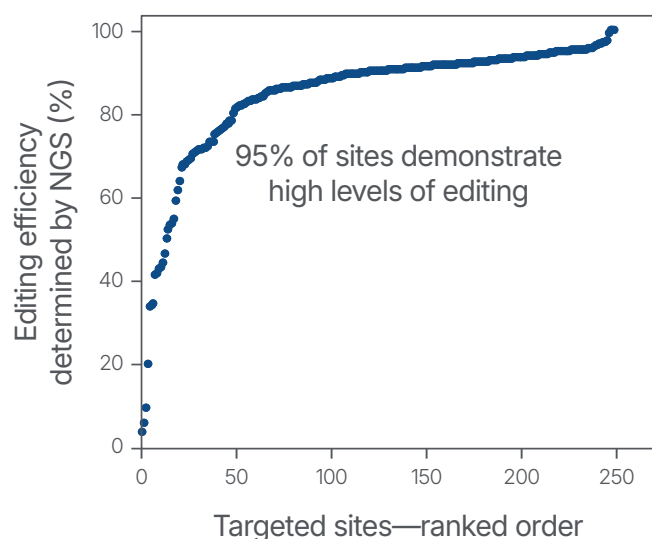



Figure 1. Alt-R CRISPR-Cas9 sgRNAs result in high editing rates in Jurkat cells. Ribonucleoprotein (RNP) complexes were formed with Alt-R S.p. WT Cas9 Nuclease V3, combined with Alt-R Cas9 sgRNAs synthesized for 255 randomly selected Cas9 guide RNA sites ($n = 1$ per site) across the human genome. RNP complexes (4 μ M) were delivered into Jurkat cells (human T lymphocyte-derived cancer cells) via a Nucleofector™ system (Lonza) in the presence of Alt-R Cas9 Electroporation Enhancer. Genome editing efficiencies were determined by target amplification followed by next generation sequencing on an Illumina® instrument.

Better QC for more accurate purity


A1. High-fidelity purity assessment

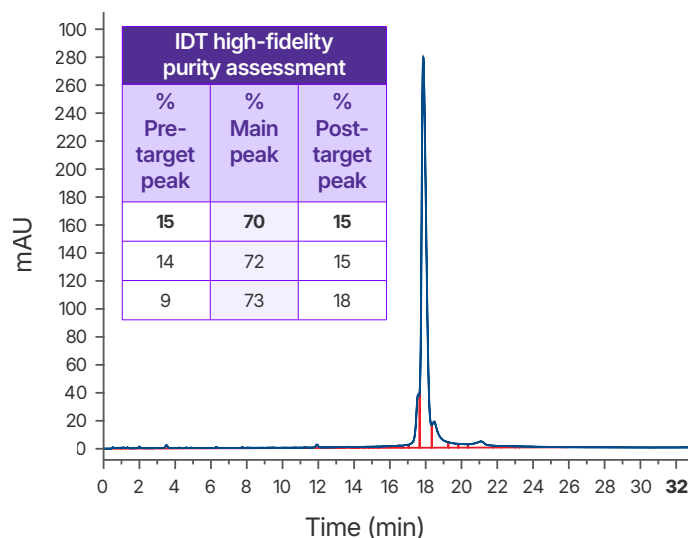

A2. Low-fidelity purity assessment

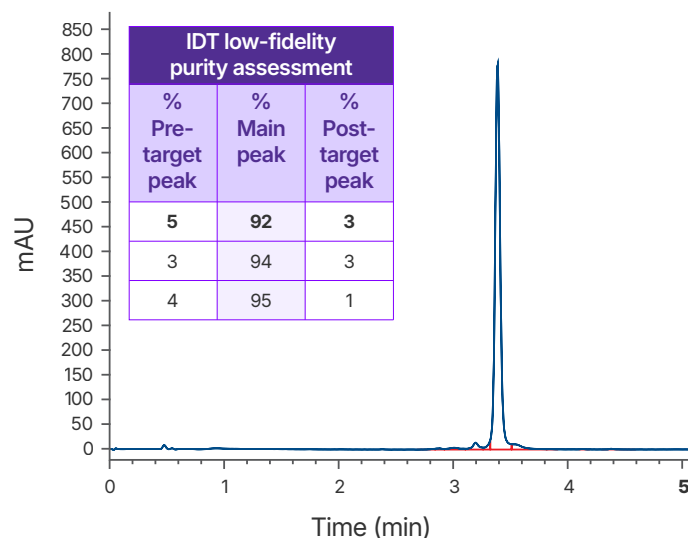

B1. Purity comparison via IDT's high-fidelity UHPLC

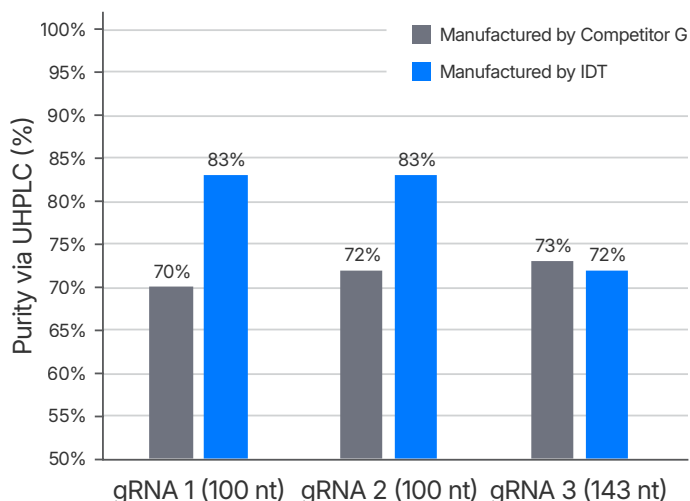

B2. Average HPLC gRNA purity comparison

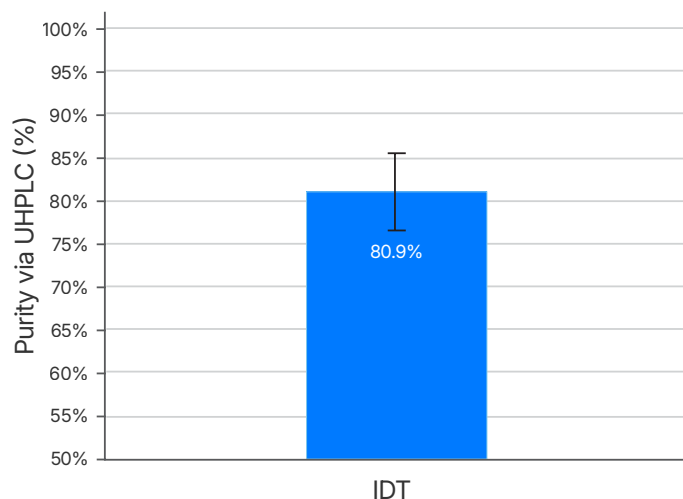


Figure 2. IDT's gRNA UHPLC purity assessment method vs a leading competitor. Left chromatogram (A1) represents purity of competitor G sgRNA using IDT's high-fidelity QC method. Right chromatogram (A2) is purity as reported by a low-fidelity QC method performed at IDT which does not allow enough time for sample separation and indicates artificially high purity. This low-fidelity method was developed to demonstrate the importance of a longer run time for purity accuracy. It was repeated for 3 guides ranging from 100–143 nt (one representative chromatogram shown). (B1) Average gRNA purity of three gRNAs of identical sequence from a competitor and IDT using IDT's **high-fidelity QC** to eliminate inflated purity estimates. (B2) Average purity of IDT HPLC gRNAs with a range of chemical modification patterns $n = 102$.

Increased stability and efficiency with chemical modifications

Available modifications*

RNA, 2'O-Methyl, phosphorothioate bonds, DNA, 2' Fluoro, Affinity Plus™ Locked Nucleic Acids

* Not all modifications are available in combination and on all types of guide RNAs.

For more information, visit [idtdna.com/CRISPR](https://www.idtdna.com/CRISPR)



For Research Use Only. Not for diagnostic procedures. Unless otherwise agreed to in writing, IDT does not intend these products to be used in clinical applications and does not warrant their fitness or suitability for any clinical diagnostic use. Purchaser is solely responsible for all decisions regarding the use of these products and any associated regulatory or legal obligations.

© 2025 Integrated DNA Technologies, Inc. All rights reserved. For specific trademark and licensing information, see www.idtdna.com/trademarks. Doc ID: RUO22-1220_002 10/25