

Resuspension of duplexed oligonucleotides (<50 nmol yield*)

resuspension protocol

A protocol for resuspending dried, annealed oligos, including Dicer-Substrate siRNAs (DsiRNAs)

1. Centrifuge tubes before opening to ensure duplexed oligos are at the bottom of the tube.
2. Resuspend duplexed oligos in Nuclease-Free Water (catalog # 11-04-02-01) to make a stock solution (concentration $\geq 100 \mu\text{M}$). For example:
3. Make further dilutions ($< 100 \mu\text{M}$) using a buffer containing 100 mM Na^+ or K^+ (e.g., a Nuclease-Free Duplex Buffer [100 mM Potassium Acetate; 30 mM HEPES, pH 7.5, catalog # 11-05-01-03]). For example:

Duplexed oligo amount	Nuclease-Free Water (100 μM final concentration)
2 nmol	20 μL
10 nmol	100 μL
25 nmol	250 μL
50 nmol	500 μL

Final concentration	100 μM duplexed oligo (from Step 2)	Buffer (containing 100 mM Na^+ or K^+)
50 μM	20 μL	20 μL
20 μM	20 μL	80 μL
10 μM	10 μL	90 μL

To calculate other dilutions, use the online IDT Dilution Calculator at <https://www.idtdna.com/Calc/Dilution/>.

Visit www.idtdna.com/protocols to verify that you are using the most current version of this protocol.

* For instructions for resuspending duplexed oligos of $\geq 50 \text{ nmol}$ yield, turn card over. Salts will be present following annealing and dry-down processes at IDT. To maintain suitable salt concentrations for the duplex structure of your product, we recommend these resuspension protocols.

For Research Use Only. Not for use in diagnostic procedures. Unless otherwise agreed to in writing, IDT does not intend these products to be used in clinical applications and does not warrant their fitness or suitability for any clinical diagnostic use. Purchaser is solely responsible for all decisions regarding the use of these products and any associated regulatory or legal obligations.

© 2023 Integrated DNA Technologies, Inc. All rights reserved. Trademarks contained herein are the property of Integrated DNA Technologies, Inc. or their respective owners. For specific trademark and licensing information, see www.idtdna.com/trademarks. Doc ID: RUO-10214-1122 01/23



WWW.IDTDNA.COM


INTEGRATED DNA TECHNOLOGIES

Resuspension of duplexed oligonucleotides (≥ 50 nmol yield*)

resuspension protocol

A protocol for resuspending dried, annealed oligos, including Dicer-Substrate siRNAs (DsiRNAs)

1. Centrifuge tubes before opening to ensure duplexed oligos are at the bottom of the tube.
2. Resuspend duplexed oligos in Nuclease-Free Water (catalog # 11-04-02-01) to make a stock solution (volume ≤ 500 μL). For example:
3. Make further dilutions (>500 μL) using a buffer containing 100 mM Na^+ or K^+ (e.g., a Nuclease-Free Duplex Buffer [100 mM Potassium Acetate; 30 mM HEPES, pH 7.5, catalog # 11-05-01-03]). For example:

To calculate other dilutions, use the online IDT Dilution Calculator at <https://www.idtdna.com/Calc/Dilution/>.

Visit www.idtdna.com/protocols to verify that you are using the most current version of this protocol.

Final concentration	Duplexed oligo amount	Nuclease-Free Water
200 μM	100 nmol	500 μL
500 μM	250 nmol	500 μL

Final concentration	200 μM duplexed oligo (from Step 2)	Buffer (containing 100 mM Na^+ or K^+)
50 μM	25 μL	75 μL
10 μM	5 μL	95 μL

Final concentration	500 μM duplexed oligo (from Step 2)	Buffer (containing 100 mM Na^+ or K^+)
50 μM	10 μL	90 μL
10 μM	2 μL	98 μL

* For instructions for resuspending duplexed oligos of <50 nmol yield, turn card over. Salts will be present following annealing and dry-down processes at IDT. To maintain suitable salt concentrations for the duplex structure of your product, we recommend these resuspension protocols.

For Research Use Only. Not for use in diagnostic procedures. Unless otherwise agreed to in writing, IDT does not intend these products to be used in clinical applications and does not warrant their fitness or suitability for any clinical diagnostic use. Purchaser is solely responsible for all decisions regarding the use of these products and any associated regulatory or legal obligations.

© 2023 Integrated DNA Technologies, Inc. All rights reserved. Trademarks contained herein are the property of Integrated DNA Technologies, Inc. or their respective owners. For specific trademark and licensing information, see www.idtdna.com/trademarks. Doc ID: RUO-10214-1122 01/23



WWW.IDTDNA.COM


INTEGRATED DNA TECHNOLOGIES